**Operator Instructions**

2 – 11

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12 – 21

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22 – 31

**Manual Del Operador**

32 – 41

**Notes**

42 – 43

# Biohit Prospenser Operator Instructions

For capacities of


2.5ml

5ml

10ml

30ml and

50ml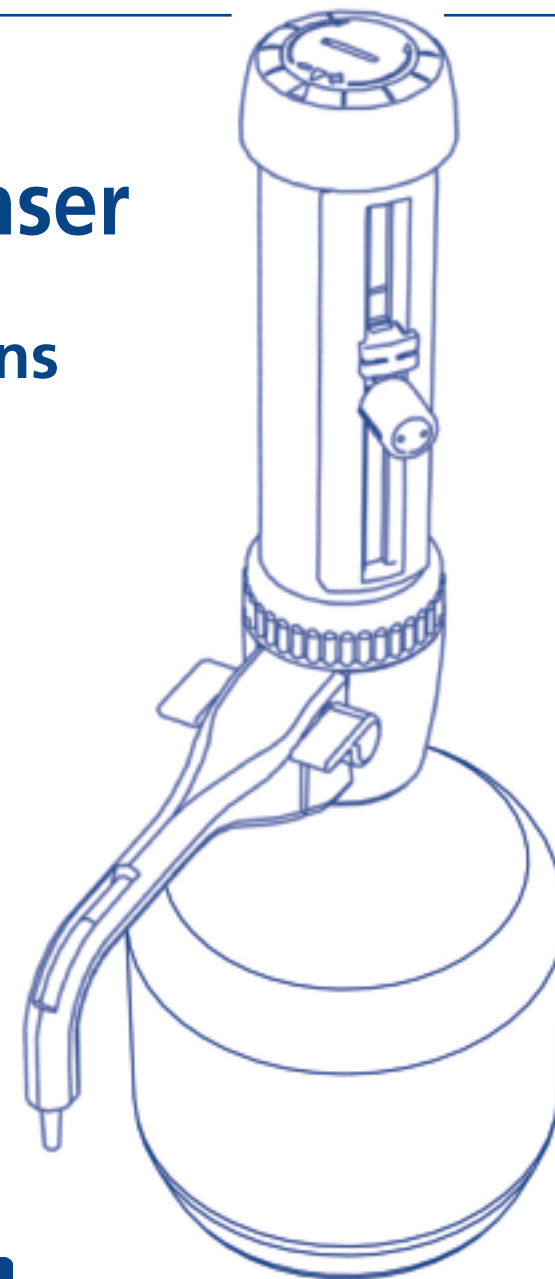


<p><b>Biohit Plc. Headquarters</b> Laippatie 1, 00880 Helsinki, Finland Tel: +358-9-773 861 Fax: +358-9-773 86200 E-mail: info@biohit.com www.biohit.com</p>	<p><b>CHINA</b> Finland Biohit Co., Ltd. Tel: +86-21-6248 5589 Fax: +86-21-6248 7786 E-mail: info.china@biohit.com</p>	<p><b>FRANCE</b> Biohit S.A. Tel: +33-1-3088 4130 Fax: +33-1-3088-4102 E-mail: commercial.france@biohit.com</p>
<p><b>GERMANY</b> Biohit Deutschland GmbH Tel: +49-6003-82 820 Fax: +49-6003-82 8222 E-mail: info@biohit.de</p>	<p><b>JAPAN</b> Biohit Japan Co., Ltd. Tel: +81-3-5822 0021 Fax: +81-3-5822 0022 E-mail: sales@biohit.co.jp</p>	<p><b>RUSSIA</b> Biohit OOO. Tel: +7-812-327 5327 Fax: +7-812-327 5323 E-mail: main@biohit.ru</p>
<p><b>U.K.</b> Biohit Ltd. Tel: +44-1803-315 900 Fax: +44-1803-315 530 E-mail: info@biohit.co.uk</p>	<p><b>U.S.A.</b> Biohit Inc. Tel: +1-732-922-4900 Fax: +1-732-922-0557 E-mail: pipet@biohit.com</p>	

## Contents

General Safety Instruction	This page
Restriction of Use	3
Before Using the Biohit Prospenser	3
Assembly	3
Operating Instructions	3
Fine Adjustment	6
Safety Precautions	6
Maintenance/Cleaning	7
Sterilisation/Autoclaving	7
Fluid-Path disassembly procedure	8
Troubleshooting	10
Accessories	11

### General Safety Instruction

- Never leave the Biohit Prospenser on your work bench with the piston Barrel full of liquid without securing the dispense anti-drip tap into 'Closed' position.
- Always check that the dispense anti-drip tap is in the 'Open' position before attempting to dispense.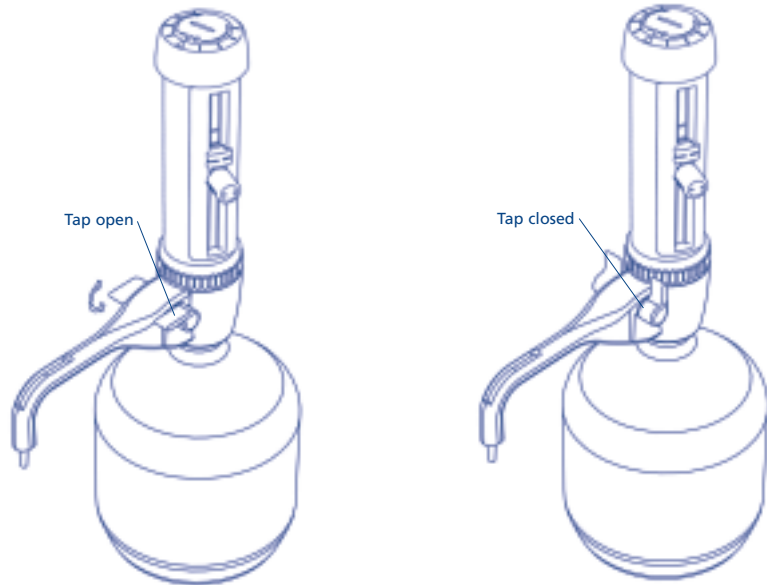

### Restriction of Use

NEVER use the Biohit Prospenser with:

- Liquids which are not compatible with PTFE, PVDF, FEP, Borosilicate glass, Alumina ceramic or might attack Platinum-Iridium
  - Hydrofluoric acid
  - Liquids which contain solid particles
- Temperature limits of Biohit Prospenser and reagent are 15°C to 40°C

### Before Using the Biohit Prospenser

Check that the instrument has not been damaged in transit.

### Assembly

The Biohit Prospenser is packed with the dispense tube attached and the inlet feed tube removed. The 300mm length of PTFE inlet tube provided should be trimmed to fit your particular reservoir. The bottom end should be cut diagonally at a length that is close to the bottom of the container when the Biohit Prospenser is assembled to the container.

The assembled Biohit Prospenser is screwed to the reservoir using gentle hand torque applied to the Biohit Prospenser's threaded platform base only. Removal should also be by means of hand torque applied to this same base.

Do not operate the piston until the unit is safely and fully mounted on the reservoir bottle.

Three adaptors are supplied to suit containers with a 33mm, 38mm or 45mm screw neck.

### Operating Instructions

#### Priming Reservoir


Fill the reservoir to approximately 50 mm below side neck aperture.

Place a container under the Biohit Prospenser dispense tube (14). Move the cursor (6) to the bottom of the scale to enable free movement of the piston. To move the cursor, simply press button (7) on the cursor sideways and slide up or down. The cursor will lock automatically into position as soon as you release the button. Prime the unit with a few gentle up and down strokes, taking the piston right down and lifting it up about 25 mm. Repeat until steady, bubble free flow is dispensed.


#### Dispensing

After priming, the piston (4) should be left in the fully down position. To dispense a precise volume of fluid, adjust the cursor (6) so that the cursor line aligns to the volume required. Draw in the required volume of fluid by lifting the piston slowly until the stop. To dispense, push the piston assembly down fully, slowly, until it stops, allowing the liquid to be collected in an appropriate vessel. When the Biohit Prospenser is not in use it is recommended that the anti-drip tap is closed.

## Items & Spares list Biohit Prospenser 2.5ml/5ml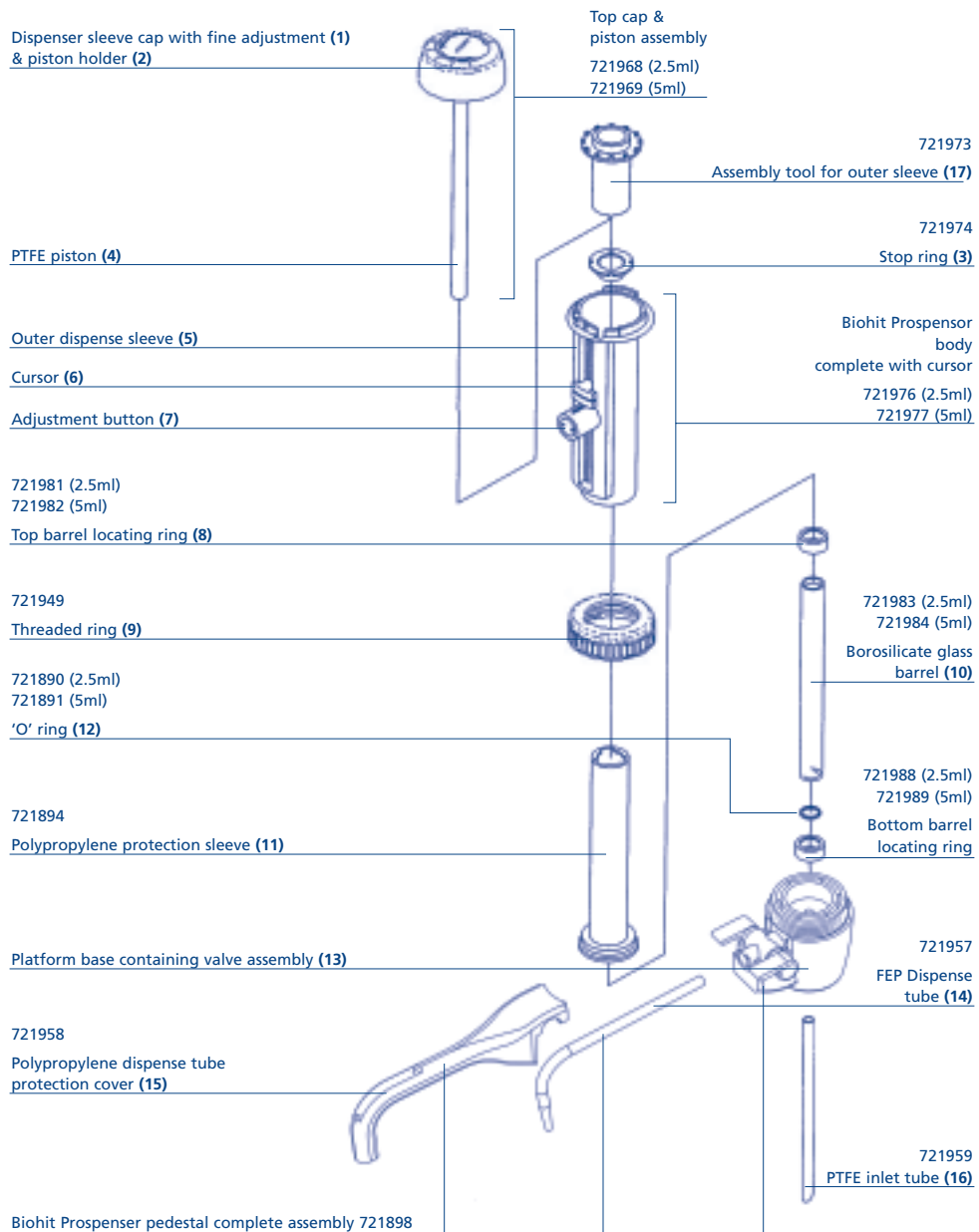

## Items & Spares list Biohit Prospenser 10ml/30ml/50ml

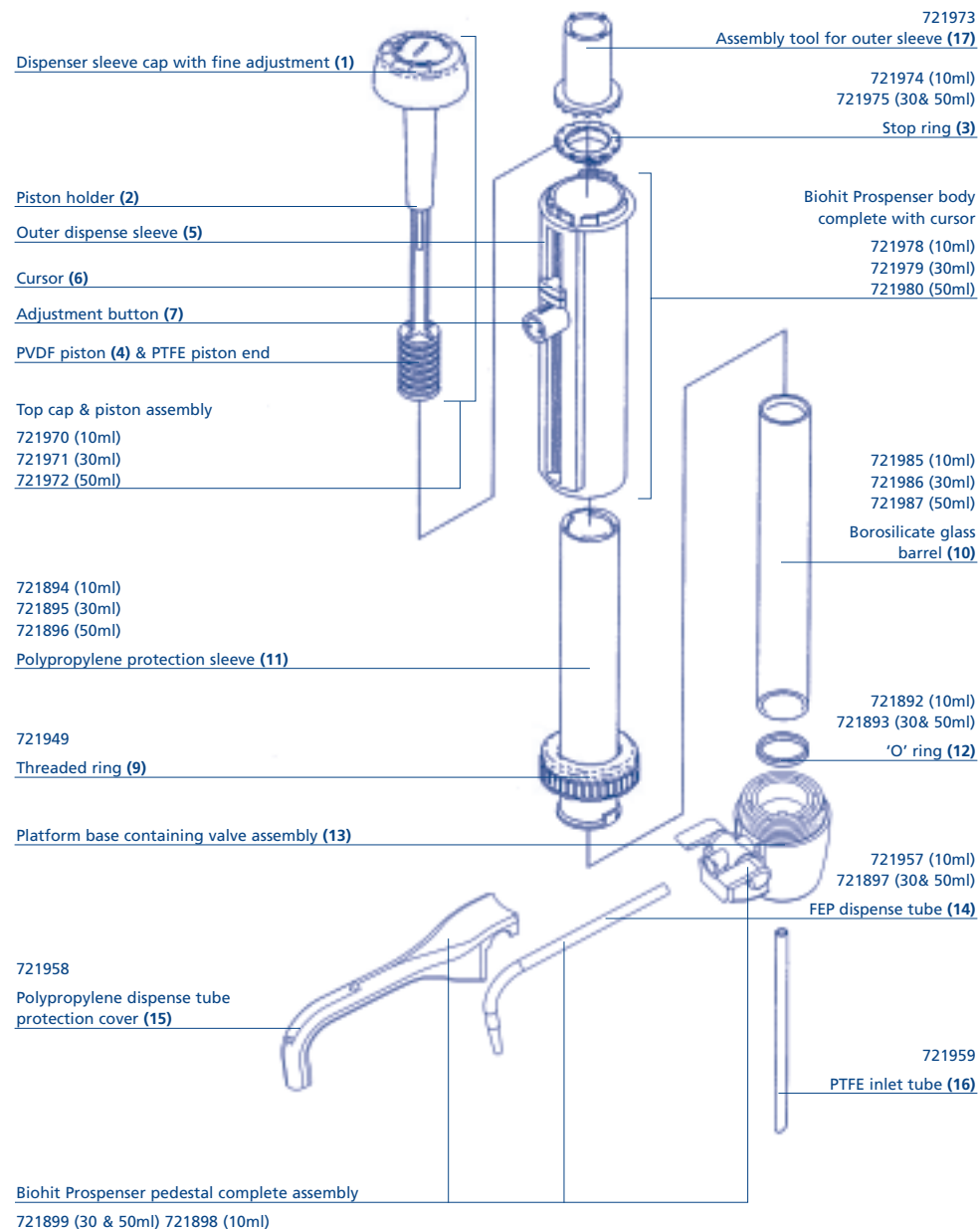
## Fine Adjustment

The figures quoted on our Test Certificate apply to the calibration of the Biohit Prospenser at full volume. The Biohit Prospenser adjustment cap enables the user to fine calibrate for accuracy. Using the adjustment disc tool that is held on the rear of the Biohit Prospenser body, change the factory calibration by turning the slot on the printed top cap insert:-

- Anti-clockwise to increase volume
- Clockwise to decrease volume

The graduations on the adjustment ring increase or decrease by the following volumes:-

Biohit Prospenser 50ml	0.1ml per line
Biohit Prospenser 30ml	0.1ml per line
Biohit Prospenser 10ml	33µl per line
Biohit Prospenser 5ml	17µl per line
Biohit Prospenser 2.5ml	8µl per line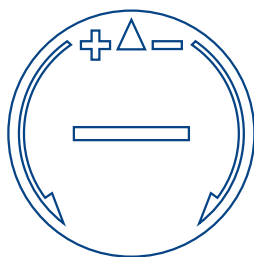

## Safety Precautions

Use the utmost caution when dispensing caustic, radioactive or hazardous chemicals.

- 1 Follow safety regulations (eg. protective clothing, goggles etc.)  
If in doubt, consult your safety officer.
- 2 Follow the operating instructions.
- 3 Only use the instrument for its proper purpose and within the limits of its materials. If in doubt, please consult the manufacturer/distributor.
- 4 Regularly inspect the instrument for leakage. Before use it is sensible to check connections.
- 5 Do not use force. Damage to parts may cause leaks.
- 6 The temperature of the instrument and reagent to be dispensed must not exceed 40°C.
- 7 Please note,
  - (i) only the manufacturer's original parts must be used.
  - (ii) materials coming into contact with the liquid to be dispensed are, borosilicate glass, PTFE, Alumina ceramic, FEP, PVDF and Platinum Iridium.

## Maintenance/Cleaning

Note; All maintenance should be carried out wearing suitable eye protection and protective clothing. If in doubt, consult your safety officer.

1. Make sure that the Biohit Prospenser is completely empty and turn the anti-drip tap to 'Open' position.
2. Place the instrument into an empty sink together with its reservoir. Unscrew the threaded platform base from the reservoir and lift the dispenser's intake tube carefully out of the reservoir, whilst tapping it against the reservoir's aperture to shake off any droplets from the intake tube.
3. Hold the dispense tube nozzle over the aperture of the reservoir and apply gentle piston strokes in order to return any syringe contents into the reservoir.
4. Flush out remaining syringe contents with distilled water or a suitable solvent. This will preserve the smooth action of the piston and free action of the inlet and outlet valves. If the inlet valve does stick and is not freed by flushing, it may be freed by gently inserting a thin rod into the inlet aperture and gently pushing the ball off its seating. Re-check unit for operation.
5. Empty the instrument completely after cleaning, and flush through with distilled water.


## Autoclaving

Steam sterilization is permissible at 121°C 2 bar after the normal cleaning procedure is carried out. The cursor should be locked in the centre of the adjustment slot before autoclaving.

Place the instrument on a cloth in order to avoid contact with metal surface. Steam sterilize the unit in its assembled status but with the threaded ring completely loosened. Steam sterilization of the piston outside the glass barrel will damage it. Cool slowly back to ambient temperature before use.

## Chemical sterilisation

The fluid path components of the Biohit Prospenser can be soaked overnight in a dilute (1:1000) solution of Sodium Hypochloride. (See page 8 for disassembly procedure).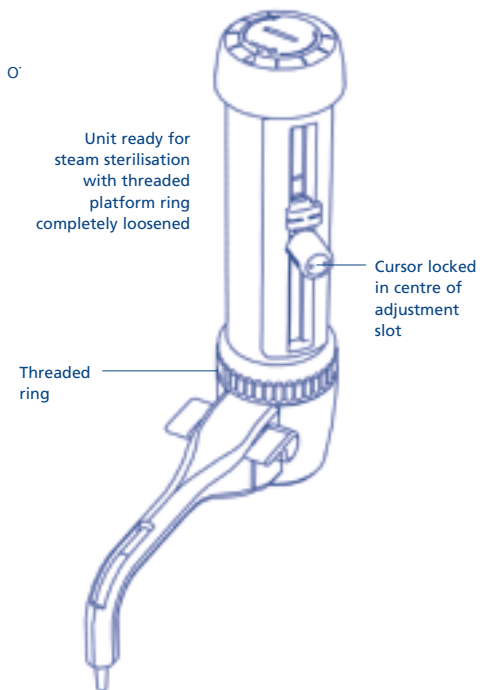


## Disassembly Options/Cleaning

### **WARNING Do not use force in assembly or disassembly.**

Disassembly should only be undertaken **after** the unit has been cleaned using the recommended cleaning procedure. Wear protective clothing and goggles during disassembly.

- 1 Move the cursor (6) down the scale and undo the dispense sleeve cap (1) by turning anti-clockwise and lifting.
- 2 The piston holder (2) is now exposed and the piston (4) is ready for removal. Remove the piston assembly carefully.
- 3 Clean the piston's surface with a suitable cleaning solution. Do not use hard tools to scrape off residue of reagent from the piston as this will damage the surface and will not seal with the glass barrel when reassembled
- 4 If it is necessary to clean or replace the graduated outer dispense sleeve (5) of your Biohit Prospenser, this is carried out with the help of the supplied assembly tool (17).
  - (a) Loosen the Polypropylene threaded ring (9) on the base by approximately one turn.
  - (b) Using the assembly tool (17) , undo the top stop ring (3) by turning clockwise and lifting it out. You can now lift off the dispense sleeve (5).
  - (c) To remove the Polypropylene protection sleeve (11) and borosilicate glass barrel (10), fully undo the threaded ring (9), remove the polypropylene protection sleeve (11) and gently lift out the glass barrel (10) for replacement or cleaning.

**Note: Bevelled inside edge of barrel must be at the top end when fitted.**
  - (d) Removal of FEP dispense tube (14); Firstly turn the anti-drip tap to the closed position. Then remove dispense tube protection cover (15) as shown on page 9 in an upward direction allowing dispense tube (14) to be pulled out of anti-drip shut-off valve assembly in direction indicated.
  - (e) Reassemble in reverse, make sure the front end of the dispense tube is clipped into the dispense tube protection cover (15) from the underside.
  - (f) Reassemble the remaining components following the above points in reverse order, making sure that the FEP 'O' ring (12) is well located into its platform recess. **Take special care not to damage the piston when inserting it into the glass barrel.**
  - (g) Before tightening the threaded ring (9) fully onto the threaded platform base (13), rotate the Polypropylene protection sleeve (11) so that the tooth on its base locates properly into the notch on the threaded ring (9).
  - (h) After reassembling the instrument, prime with distilled water to ensure that assembly has been correctly followed and the piston is working smoothly. Check that no leaks occur.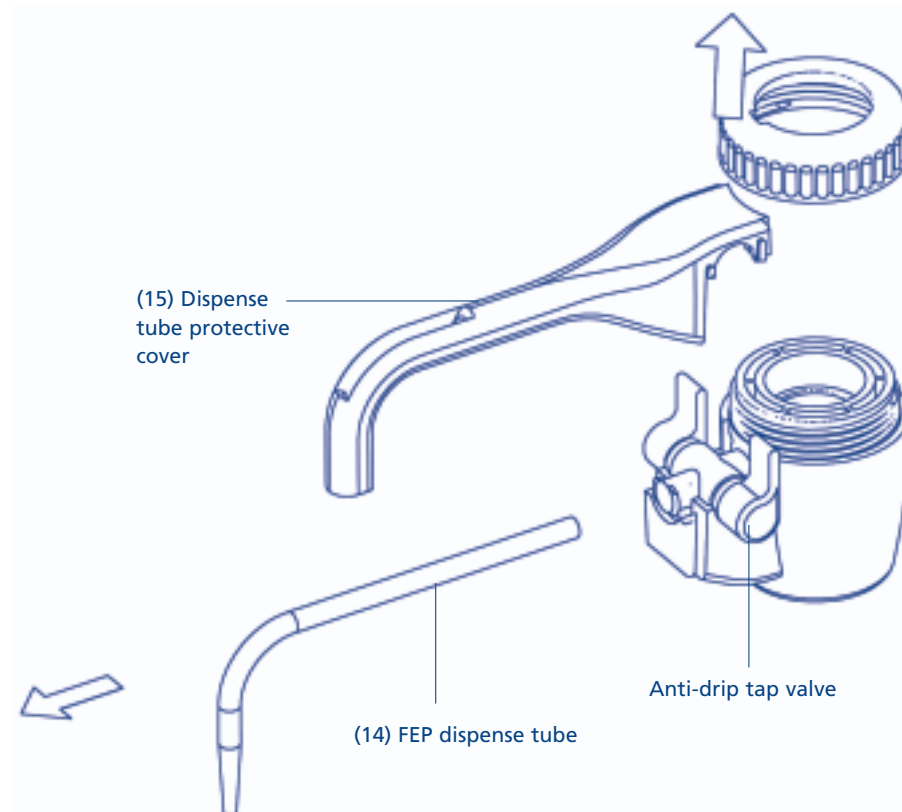


Problem	Possible Cause	Remedy
Air bubbles appear in discharge tube	Liquid reservoir is empty Too fast filling action Glass barrel is not sealing against FEP 'O' ring Leaking piston Leaking discharge valve	Refill reservoir and prime unit Fill and dispense more slowly Unscrew the threaded platform ring and make sure that the FEP 'O' ring is properly fitted into its recess Clean piston. If problem persists, replace piston Clean unit by flushing through - if problem persists, replace platform base
Barrel does not fill with liquid	Inlet tube not fitted correctly Inlet valve stuck Glass barrel is not sealing against the FEP 'O' ring	Connect inlet tube correctly Free inlet valve by inserting a thin rod into the inlet aperture Unscrew the threaded ring and make sure that the FEP 'O' ring is properly fitted into its recess
Dispensing not possible	Blocked dispense tube Discharge valve stuck	Disassemble the dispense tube and flush through with cleaning fluid Clean unit by immersing platform in cleaning fluid (taking care to protect LCD) - if problem persists, replace platform base
Wrong dispense volume	Instrument not calibrated Leaking valves	See page 6 'User Calibration' Clean platform base - if problem persists, replace
Liquid appears between glass barrel and Polypropylene protection sleeve	Glass barrel is not sealing against the FEP 'O' ring Damaged FEP 'O' ring	Unscrew the threaded ring and make sure that the FEP 'O' ring is properly fitted into its recess Replace FEP 'O' ring

## Accessories

### Adaptors

A range of five adaptors for fitting your dispenser to the reservoir are available.

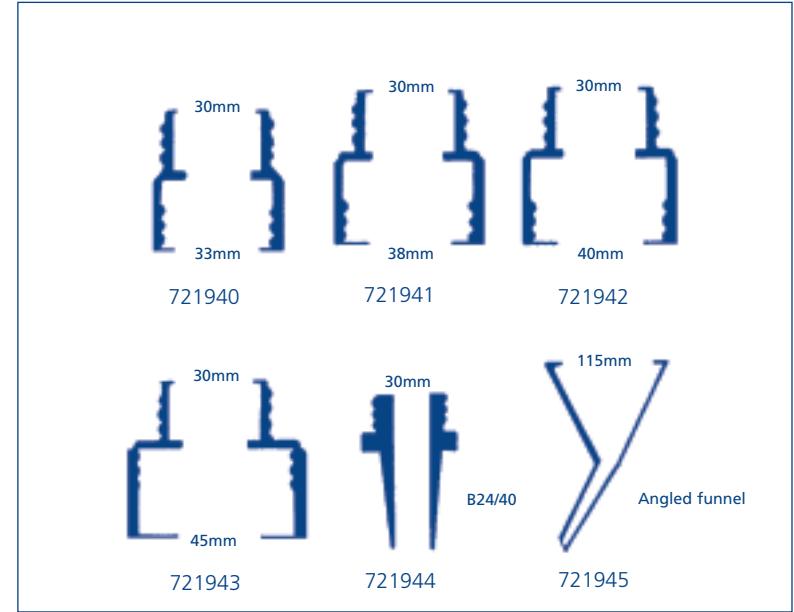### Specification

Accuracy  $\pm 0.3\%$  on maximum delivery and a precision  $\leq 0.1\% CV$  using distilled water at 20°C.

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